|                        |                          |
|------------------------|--------------------------|
| <b>PRECISTERM 2L.</b>  | <b>6000137 / 6000385</b> |
| <b>PRECISTERM 5L.</b>  | <b>6000138 / 6000386</b> |
| <b>PRECISTERM 12L.</b> | <b>6000140 / 6000387</b> |
| <b>PRECISTERM 20L.</b> | <b>6000141 / 6000388</b> |
| <b>PRECISTERM 45L.</b> | <b>6000142 / 6000389</b> |

**BAÑOS TERMOSTÁTICOS PARA AGUA Y ACEITE**  
**THERMOSTATIC BATHS FOR WATER AND OIL**

## INDICE

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## GENERAL INFORMATION

1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.

2) Do not install or use the equipment without reading this handbook before.

3) This handbook must always be attached to the equipment and it must be available for all users.

4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.

**5) IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**

6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 2009/104/CE, the manufacturer is not responsible for the damage that can occur.

7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

## PACKING LIST

The standard equipment consist of the following components:

| Description           | Code    | Quantity |
|-----------------------|---------|----------|
| Precistern 2L.        | 6000137 | 1        |
| Precistern 2L. 200°C  | 6000385 | 1        |
| Precistern 5L.        | 6000138 | 1        |
| Precistern 5L. 200°C  | 6000386 | 1        |
| Precistern 12L.       | 6000140 | 1        |
| Precistern 12L. 200°C | 6000387 | 1        |
| Precistern 20L.       | 6000141 | 1        |
| Precistern 20L. 200°C | 6000388 | 1        |
| Precistern 45L.       | 6000142 | 1        |
| Precistern 45L. 200°C | 6000389 | 1        |
| Instruction manual    | 80003   | 1        |

## ACCESSORIES

(The following components, at user's request, must be fitted in factory)

| Denomination | Code |
|--------------|------|
|--------------|------|

|                                                                   |         |
|-------------------------------------------------------------------|---------|
| REMOVABLE LEVEL<br>For baths to be used without perforated cover. | 6001400 |
|-------------------------------------------------------------------|---------|

|                                              |         |
|----------------------------------------------|---------|
| FIXED LEVEL<br>It must be fitted in factory. | 6000123 |
|----------------------------------------------|---------|

|                                               |         |
|-----------------------------------------------|---------|
| THERMOMETER<br>To adapt with bosshead (120°C) | 1000120 |
| To adapt with bosshead (200°C)                | 1000200 |

|                                                    |         |
|----------------------------------------------------|---------|
| THERMOMETER<br>To be fitted at gable cover (120°C) | 1000872 |
| To be fitted at gable cover (200°C)                | 1000873 |

|                                  |         |
|----------------------------------|---------|
| SUPPORT BOSSHEAD FOR THERMOMETER | 6000896 |
|----------------------------------|---------|

Set opening up to 40mm.  
In AISI 304 stainless steel.

| For bath code                                                                                                                                                                      |          | Cover with holes and reducing discs 105, 80, 60 and 37 mm diam, with a hole for the thermometer in AISI 304 stainless steel | Gable cover in AISI 304 stainless steel | Perforated shelf in AISI 304 stainless steel |
|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|----------|-----------------------------------------------------------------------------------------------------------------------------|-----------------------------------------|----------------------------------------------|
| 6000137                                                                                                                                                                            | 6000385  | 6003231 (1 hole)                                                                                                            | No                                      | 6000011                                      |
| 6000138                                                                                                                                                                            | 6000386  | 6003232 (2 holes)                                                                                                           | 6000710                                 | 6000022                                      |
| 6000140                                                                                                                                                                            | 6000387  | 6003234 (4 holes)                                                                                                           | 6000711                                 | 6000044                                      |
| 6000141                                                                                                                                                                            | 6000388  | 6003236 (6 holes)                                                                                                           | 6000712                                 | 6000066                                      |
| 6000142*                                                                                                                                                                           | 6000389* | 2x6003236 (6 holes) = (12 holes)                                                                                            | 6000714                                 | 6000088                                      |
| *Because of the large surface of baths code 6000142 and 6000389, to work at temperatures above 50°C, the use of cover with holes and reducing discs or gable cover is recommended. |          |                                                                                                                             |                                         |                                              |

## TECHNICAL FEATURES

**Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.**

| Code     | Maximum temperature °C | Capacity litres | Tank dimensions cm |        |       | Overall dimensions cm |        |       | Consumption W | Weight Kg |
|----------|------------------------|-----------------|--------------------|--------|-------|-----------------------|--------|-------|---------------|-----------|
|          |                        |                 | Depth              | Length | Width | Depth                 | Length | Width |               |           |
| 6000137  | 110                    | 2               | 15                 | 15     | 13    | 20                    | 20     | 27    | 390           | 3.5       |
| 6000385  | 200                    |                 |                    |        |       |                       |        |       | 640           |           |
| 6000138  | 110                    | 5               | 15                 | 30     | 14    | 20                    | 37     | 27    | 600           | 6         |
| 6000386  | 200                    |                 |                    |        |       |                       |        |       | 1110          |           |
| 6000140  | 110                    | 12              | 15                 | 30     | 30    | 20                    | 40     | 42    | 1110          | 8         |
| 6000387  | 200                    |                 |                    |        |       |                       |        |       | 1400          |           |
| 6000141  | 110                    | 20              | 15                 | 48     | 30    | 20                    | 58     | 42    | 1500          | 10        |
| 6000388  | 200                    |                 |                    |        |       |                       |        |       | 2100          |           |
| 6000142* | 110                    | 45              | 15                 | 62     | 50    | 22                    | 69     | 62    | 2400          | 18        |
| 6000389* | 200                    |                 |                    |        |       |                       |        |       | 3800          |           |


\* With drain tap.

The temperature stability is  $\pm 1$  K for baths of 100°C and  $\pm 2$  K for baths of 200°C.

## EQUIPMENT DESCRIPTION

**PRECISTERM** baths with adjustable liquid temperature from ambient +5°C up to 110°C (stability of  $\pm 1^\circ\text{C}$ ) or up to 200°C (stability of  $\pm 2^\circ\text{C}$ ), depending on the model. To use the baths, fill the tank with liquid until the level covers at least the heater element. The liquid or mixture to be employed must be adequate to the model of bath and the working temperature. As a guidance, here are some liquids:  
(To use these liquids, follow the manufacturer's instructions).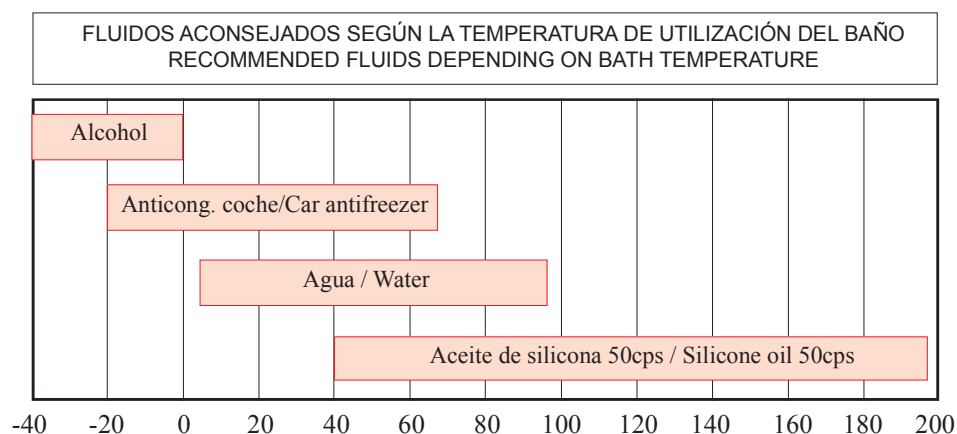

## INSTALLATION

Place the bath on a flat, horizontal and level surface. Bear in mind the weight of the liquid when choosing the support. Provide for bath drainage (bath 6000142 and 6000389).

**Be sure that the voltage supply is the same as the one indicated on the characteristics plate.**

**Do not use the equipment if it is not earthed.**

**If you change the plug bear in mind the following:**

**Blue cable: Neutral.**

**Brown cable: Phase.**

**Yellow/green cable: Earth.**

**It is recommended the use of decalcified water. DO NOT USE DEIONIZED OR DEMINERALIZED WATER as it has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of  $\text{CO}_3\text{Na}_2$  (Sodium Carbonate) per litre of water used in the bath.**


**CAUTION!!! IMPORTANT FOR YOUR SAFETY****OPERATION****CONTROL PANEL:**

1. Main switch with indicator lamp.
2. Temperature control knob.
3. Heater operation indicator lamp (amber).
4. Safety thermostat operation indicator lamp (red).

**REAR PART**

5. Safety thermostat control screw.
6. Safety thermostat manual reset button (grey).

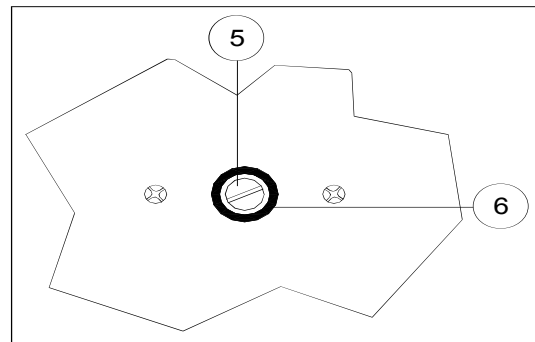**STARTING UP OPERATION :**

1. Fill the bath with liquid. Leave at least 3 cm without filling to avoid splashes.
2. Connect the bath to the mains.
3. Turn on the main switch (1). The green indicator lamp will light up.
4. Turn on the temperature control thermostat (2) from low to high according to the working temperature. The heater operation indicator lamp (amber) will light up (3).
5. After a certain time, the heater operation indicator lamp (3) will begin to operate intermittently. When the cycle of these intermittences becomes uniform, that will indicate that temperature of liquid is stabilized.

6. The thermostat scale graduation is only useful as a reference, because the real temperature is the one shown by a thermometer introduced into the liquid.

7. Be careful with liquid evaporation. The heating element must ALWAYS be covered with liquid. Replenish when necessary.

#### ADJUSTMENT FOR THE SAFETY THERMOSTAT:

The low limit of working temperature for the safety thermostat is nearly 50°C. The desired temperature of the safety thermostat must be at least 5°C above the working temperature. The adjusting process of the safety thermostat is long. For that reason, the safety thermostat is useful when the users work habitually at the same temperature.

1. Turn the safety thermostat screw (5) clockwise to maximum.
2. Select with the temperature control knob (2) the desired temperature of the safety thermostat.
3. Turn on the bath and leave it to stabilize.
4. When the bath is stabilized, slowly turn the safety thermostat screw (5) anticlockwise until it clicks and the red indicator lamp (4) lights up.
5. Wait until the bath cools to reset the safety thermostat with (6) button.
6. Select the working temperature of the bath with temperature control knob (2).

## SPARE PARTS

**To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, s.a.u.**

| DESCRIPTION                           | CODE          |         |         |         |         |         |         |         |         |         |
|---------------------------------------|---------------|---------|---------|---------|---------|---------|---------|---------|---------|---------|
| Model                                 | 6000137       | 6000385 | 6000138 | 6000386 | 6000140 | 6000387 | 6000141 | 6000388 | 6000142 | 6000389 |
| 1. Heating element                    | 39063         | 39049   | 39061   | 39050   | 39051   | 39060   | 39012   | 39059   | 39123   | 39124   |
| 2. Safety thermostat                  | 43016         | 43017   | 43016   | 43017   | 43016   | 43017   | 43016   | 43017   | 43016   | 43017   |
| 3. Control thermostat                 | 43006         | 43005   | 43006   | 43005   | 43006   | 43005   | 43006   | 43005   | 43006   | 43005   |
| 4. Thermostat racord + thermostat nut | 34141 + 34142 |         |         |         |         |         |         |         |         |         |
| 5. Mains cable                        | 7024          |         |         |         |         |         |         |         |         |         |
| 6. Omron contactor                    | No            |         |         |         |         |         |         |         |         | 13012   |

## MAINTENANCE

**Before removing the cover, disconnect the equipment from the mains.**

**The control element and the heating element must only be manipulated by authorized personnel.**

**The distilled or demineralized water has a great tendency to rust. To prolong the life of the bath and avoid rusting of the parts, dissolve 1 gr of  $\text{CO}_3\text{Na}_2$  (Carbonate of Soda) per litre of water used in the bath.**

### CLEANING:

If normal tapwater is used calcareous deposits will appear in the bath and in the heating element. A heater covered with lime should not be used as it can lose heating power and deteriorate more quickly. Periodical cleaning of the inside of the tank and the heating element with a solution of 10% of 28% Chlorhydric acid or a solution of water and vinegar (Acetic acid) is advised. Rinse with clean water afterwards.

For the cleaning of the different parts use the following products:


Cleaning of stainless steel: alcohol

Cleaning of plastic: Alcohol with cotton duster.

## GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

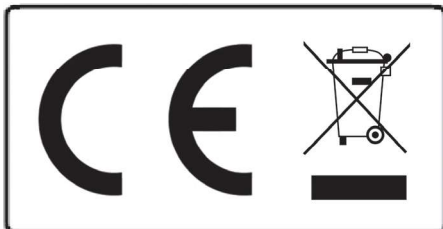### Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.