LAB BOTTLES AND ACCESORIES

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LAB BOTTLES AND ACCESORIES

Bottles for reagents are made of borosilicate glass 3.3 SIMAX. For the photosensitive reagents, the brown colored bottles are intended.

Plastic screw caps (blue) are intended for the operational temperature up to 140°C. In case of higher temperature (maximum 200°C), the red screw caps should be used.

In case of reagents, which could damage the plastic screw caps, the bottles with precise ground bottle neck and glass stopper are intended.

For storage of bigger amounts as well as for bleeding of liquids, the irrigators are delivered.

The bottles of thick-wall type are not intended for reduced or increased pressure.

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REAGENT BOTTLE WITH GL80

with blue PP cap
with blue PP outlet ring

Provided with well-legible scale and large field for description (indication of the content). The outlet ring enables for clean work (no drops). The great diameter of bottle neck makes the handling with powders or viscous substances easy.

Typical use: storage and transfer of substances.

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TP
1270

REAGENT BOTTLES WITH GL80

with blue PP cap
with blue PP pouring ring

Provided with well-legible scale and large field for description (indication of the content). The pouring ring enables for clean work (no drops). The great diameter of bottle neck makes the handling with powders or viscous substances easy. Brown colored glass provides for protection of photosensitive substances.

Typical use: storage and transfer of photosensitive substances.

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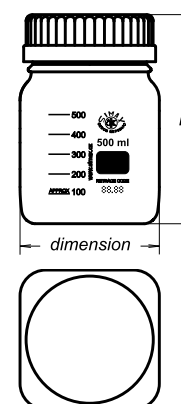
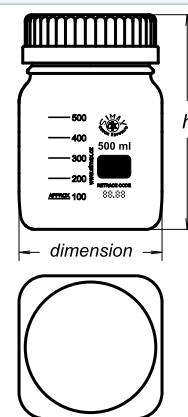
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
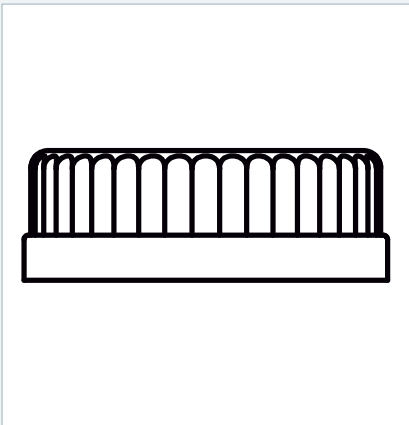
TP
1270

SCREW CAP GL 80

Plastic cap (blue) GL 80.
Maximum operational temperature
+140°C.

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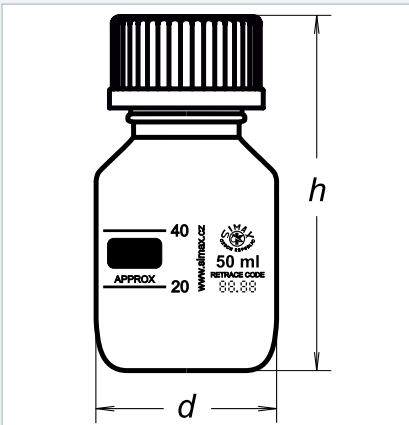
REAGENT BOTTLE

brown, with blue PP cap and blue PP outlet ring

Provided with well-legible scale and large field for description (indication of the content). The outlet ring enables for clean work (no drops). Typically used for storage and transfer of photosensitive substances.

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ISO
4796-1

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REAGENT BOTTLE WITH GL45

with red PBT cap
with red ETFE outlet ring

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
Cat. No. 2070 R

Catalog No. Katalog-Nr.	Capacity (ml) Kapazität	GL	d (mm)	h (mm)	Packing (pcs) Packung (Stück)
1632414321101	100	45	56	105	10
1632414321251	250	45	70	143	10
1632414321501	500	45	86	181	10
1632414321941	1000	45	101	230	10
1632414321951	2000	45	136	265	10
1632414321955	3800	45	160	320	4
1632414321959	5000	45	181	336	6
1632414321967	10000	45	227	416	1
1632414321969	20000	45	288	512	1

TP
1270

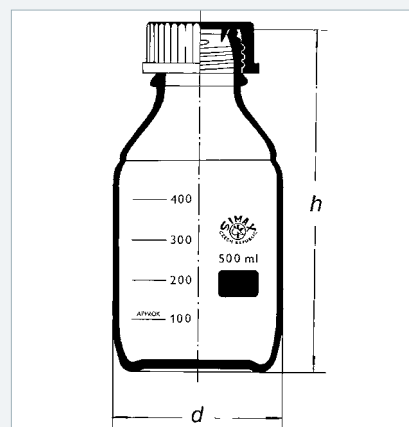ISO
4796-1

SCREW CAP

PP Maximum operating temperature
+140°C.

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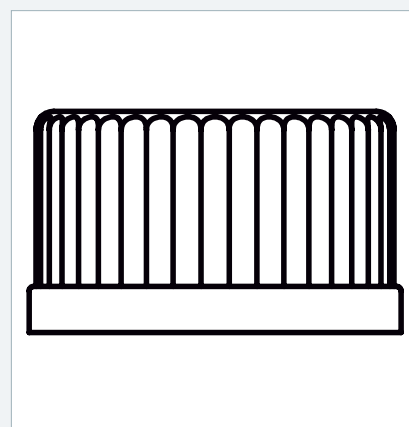
NEW
NEW
NEW

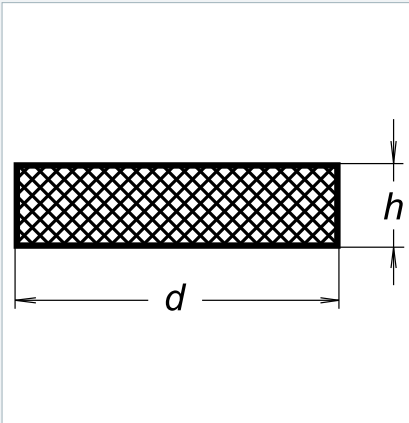
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
Note:

The red screw cap is made of PBT and maximum operating temperature is +200 °C.

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Note: The red outlet ring is made of ETFE and maximum operating temperature is +200 °C.

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SCREW FITTING

In practice, it is often necessary to interconnect piping (made of PTFE, PFA, FEP, ...) or pipes (glass, metal, plastic) with the instruments provided with GL (glass) thread. In such cases, the screw fitting presents an ideal solution.

Every screw fitting consists of the GL threaded sleeve nut, three internal parts of V-ring, conical ring, and sealing ring.

The assembling is very easy and may be performed by hand: slip the internal parts over the pipe; then insert the pipe end into the mating part (connecting port of the instrument, bottle neck, ...) and tighten the nut. In this way the nut will press both the sealing and conical ring against the mating part. At the same time the V-ring will be compressed and an absolutely tight connection is obtained, suitable even for vacuum.

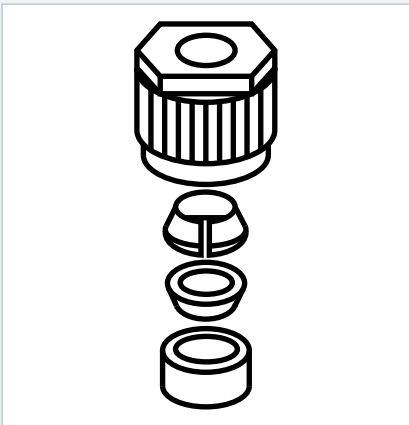
At the room temperature, the lab gaskets GL14, GL18 and GL25 bear up the pressure of maximum 10 bar.

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SCREW FITTING

Red sleeve nut (glass fibre reinforced ETFE), internal parts V-ring (ETFE), conical nut (PTFE) and sealing nut (PTFE), O-ring (for piping with diameter less than 3mm, not in contact with medium). Very good chemical resistance, suitable for temperature range -50°C to +150°C.

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SCREW FITTING

Screw fittings and fittings with GL thread for hard piping or pipes made of glass, metal, or plastic). For attachment of probes, thermometers, or cables in the reaction vessels. Ideal for the aggressive environments (aggressive vapors).

Green sleeve nut (PP) for the bottle GL45 thread and PP body. Maximum external diameter of the inserted pipe is 8,5mm. Limited chemical resistance, for operational temperatures up to +110°C. Suitable for inflow/outflow of liquids, leading-in the piping, pipes or probes into the vessels.

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SCREW CAP

with two hose connections

Green screw cap (PP) for the bottle GL45 thread plus body with hose connection (PP). The upper hose connections are bent, the lower ones are straight. Limited chemical resistance, for operational temperatures up to +110°C.

Used for inflow/outflow of liquids. Suitable for hoses (Vitron, Tygon, silicone).

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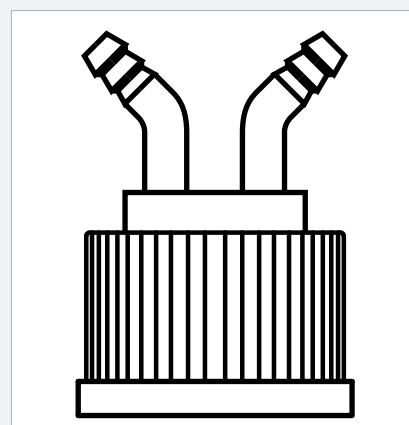
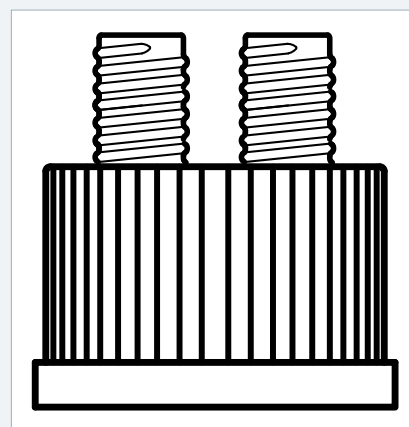
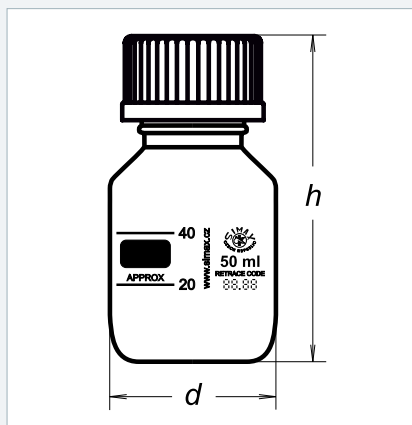
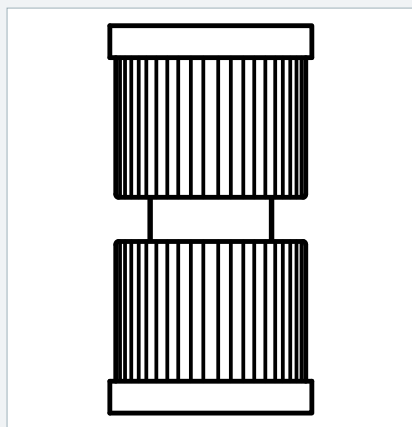
SCREW-TYPE JOINT

Two screw-type caps (PBTP, +180°C) with GL thread and PTFE/FPM gasket. Joint made of PTFE. The medium is only in contact with PTFE.

Suitable for connection of two GL threads with identical diameter.

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REAGENT BOTTLE, PLASTIC COATED

PU-polyurethane
Provided with well-legible scale and large field for description (indication of the content).

Operational temperature range (for plastic coating) -30°C to +135°C. Autoclaving: Steam sterilisation at +121°C or +134°C. The coating provides protection against scratching and rupture – ideal for transport and storage of dangerous substances and provides also protection against the light (wavelength up to approx. 380 nm). High micro-wave transmittance. Suitable for storage, transport, and safe handling with dangerous or valuable substances.

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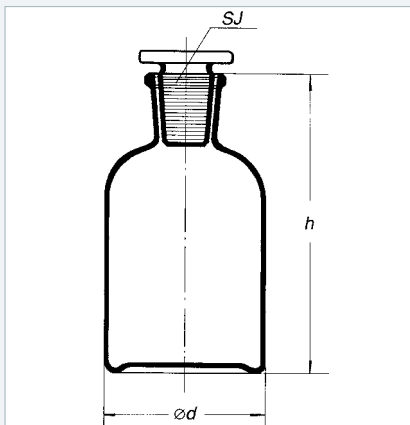
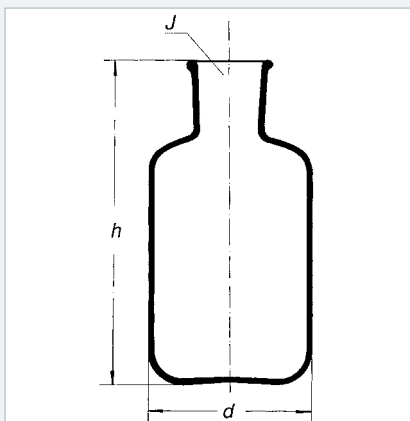
BOTTLE RESERVOIR

for injection serum

Without ground neck and stopper.
Typical use: storage of liquids (incl. the aggressive ones).

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BOTTLE REAGENT, NARROW MOUTH

with precise ground neck and glass stopper

Typical use: storage of liquids (incl. the aggressive ones).

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BOTTLE REAGENT, WIDE MOUTH

for injection serum

With precise ground neck and glass stopper.

Typical use: storage of powdery or granulated substances.

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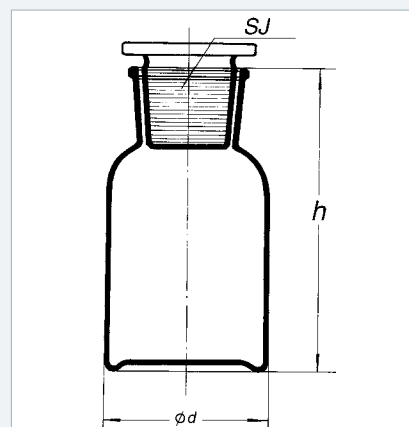
PN
70 0040

IRRIGATOR

barrel shaped, tubulated near bottom, with knob

Provided with standardized ground neck near bottom.

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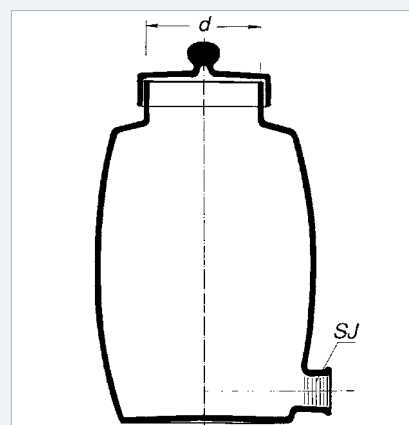
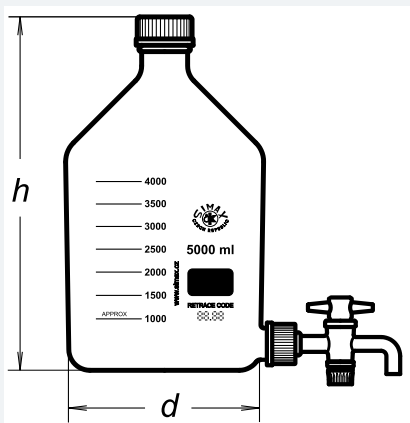
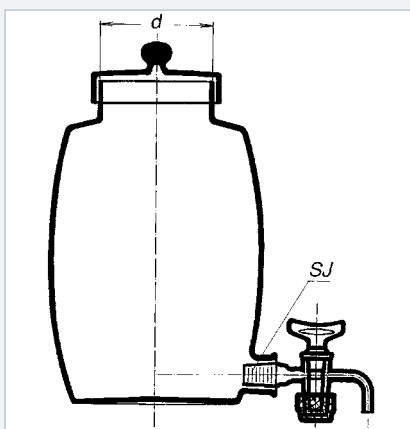
PN
70 0531

IRRIGATOR

barrel shaped, tubulated near bottom

with loose drop – on cover with knob
Provided with standardized ground neck
near bottom and bleeding cock.

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REAGENT BOTTLE

outlet GL32 + cock (PTFE)

Reagent bottle with the outlet GL32
and PTFE cock, suitable as reservoir for
liquids.

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REAGENT BOTTLE

for common use

Wide mouth - suitable for handling of powders and viscous liquids.
Typical use: storage and transport of substances.

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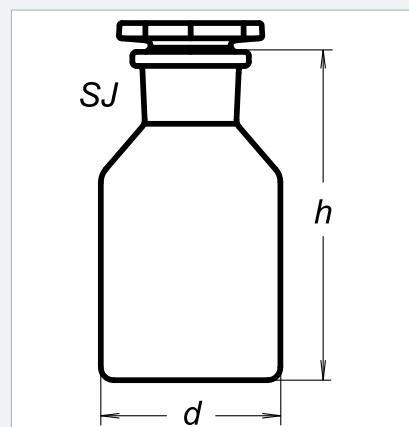
REAGENT BOTTLE

for common use, wide mouth, brown

Brown color provides for protection of photosensitive substances.
Wide mouth suitable for handling of powders and viscous liquids.
Typical use: storage and transport of photosensitive substances.

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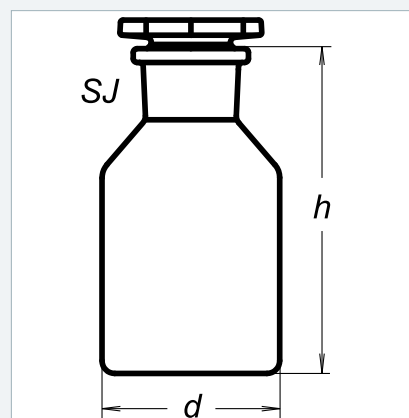
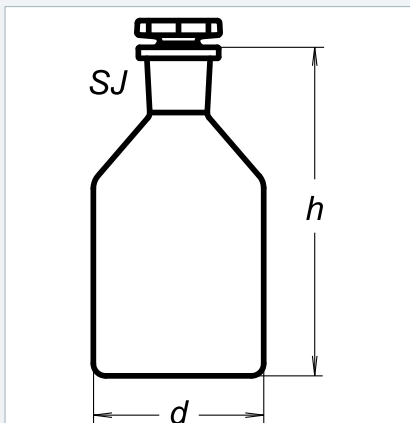
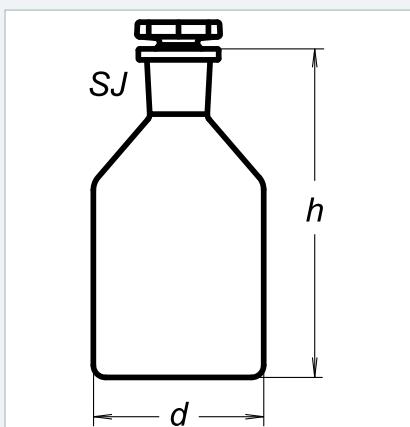
REAGENT BOTTLE

for common use

Narrow mouth, clear.
Typical use: storage and transport of liquids.

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REAGENT BOTTLE

for common use

Narrow mouth, brown.
Brown color provides for protection of photosensitive substances.
Typical use: storage and transport of photosensitive substances.

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DOSING DEVICE

Used in "in-vitro" diagnostic labs, for safe and exact dosage of agents and chemicals from the bottles.

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